CLEANJET

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Automatic nozzle for disinfecting gloved hands - AM101

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Warnings:

Read this Instruction Manual carefully before using the CLEANjet device. The use of the CLEANjet device is safe if it is carried out in compliance with the instructions contained in this Instruction Manual. This manual is the property of A.M. Instruments srl, reproduction and distribution is forbidden unless previously authorized. Manufacturer: A.M. Instruments srl - via Isonzo, 1 / C - 20812 Limbiate - MB - Italy

SUMMARY

1. Packaging content
2. Product description
3. Limits of use, residual risks and precautions
4. Installation
5. Procedures for use
6. Replacement parts
7. Warranty

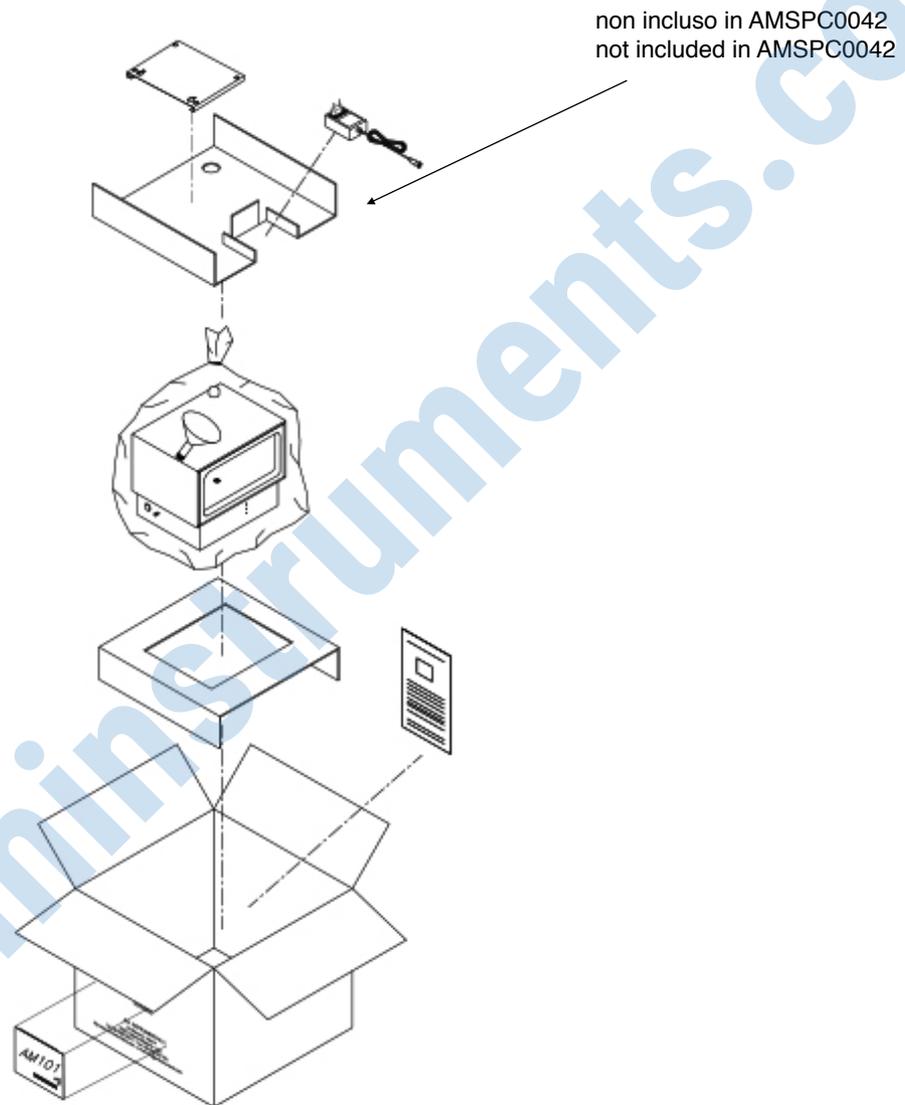
AM Instruments srl Via Isonzo, 1/C 20812 Limbiate (MB) Italy	T. +39 02 872892.1 F. +39 02 872892.00 www.aminstruments.com info@aminstruments.com	AM Instruments Tuscany Business Resort Strada di Cerchiaia 28/A 53100 Siena - Italy T. +39 02 872892.1	Sede legale Via Manzoni, 144/E-F 20811 Cesano Maderno (MB) Italy	COMPANY WITH QUALITY SYSTEM CERTIFIED BY DNV GL = ISO 9001:2015 =
--	--	--	--	---

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2. PRODUCT DESCRIPTION

2.1 – The device

CLEANjet is an automatic atomizer for the disinfection of gloved hands to be used with liquids based on alcohol. CLEANjet is not a medical device.

2.2 Technical data

3. LIMITS OF EMPLOYMENT, RESIDUAL RISKS AND PRECAUTIONS

- CLEANjet is to be used exclusively for the disinfection of gloved hands.
- Absolutely avoid introducing any other part of the body.
- Absolutely avoid introducing any object.
- Do not place the CLEANjet close to devices that emit infrared rays as they may interfere with the optical reader (7) causing the CLEANjet to operate continuously even without the introduction of gloved hands.
- Use only alcohol-based cleaning solutions without residue (eg IPA), non-foaming.
- Always ask the cleaning solution supplier for the safety data sheet that must always be available near the CLEANjet
- Properly educate all the CLEANjet users.

AM Instruments srl Via Isonzo, 1/C 20812 Limbiate (MB) Italy	T. +39 02 872892.1 F. +39 02 872892.00 www.aminstruments.com info@aminstruments.com	AM Instruments Tuscany Business Resort Strada di Cerchiaia 28/A 53100 Siena - Italy T. +39 02 872892.1	Sede legale Via Manzoni, 144/E-F 20811 Cesano Maderno (MB) Italy	COMPANY WITH QUALITY SYSTEM CERTIFIED BY DNV GL = ISO 9001:2015 =
--	--	--	--	---

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4. INSTALLATION

Installation:

The CLEANjet unit can be placed on a surface or hung on a wall.

Make sure that an electric plug compatible with that supplied with the CLEANjet is available near it. (2.2 Technical data)

If it should be hung on a wall, proceed in the following way:

- drill the wall and insert the dowels (not supplied) into the holes;
- screw the device (C) to the wall;
- hook the CLEANjet on the wall;
- take the transformer (B) being sure that the selector is in position 12V and insert the round connector (8) into the CLEANjet and connect the plug to the main socket 230V, the green light ON (1) will light up;
- the unit is now ready for the other operations.


5. INSTRUCTIONS FOR USE

The green light ON (1) placed on the front panel light up. At the same time the liquid level red light MIN (3) will also light up thus indicating the absence of cleaning solution.

If a hand is introduced a sound signal will be set off.

AM Instruments srl Via Isonzo, 1/C 20812 Limbiate (MB) Italy	T. +39 02 872892.1 F. +39 02 872892.00 www.aminstruments.com info@aminstruments.com	AM Instruments Tuscany Business Resort Strada di Cerchiaia 28/A 53100 Siena - Italy T. +39 02 872892.1	Sede legale Via Manzoni, 144/E-F 20811 Cesano Maderno (MB) Italy	COMPANY WITH QUALITY SYSTEM CERTIFIED BY DNV GL = ISO 9001:2015 =
--	--	--	--	---

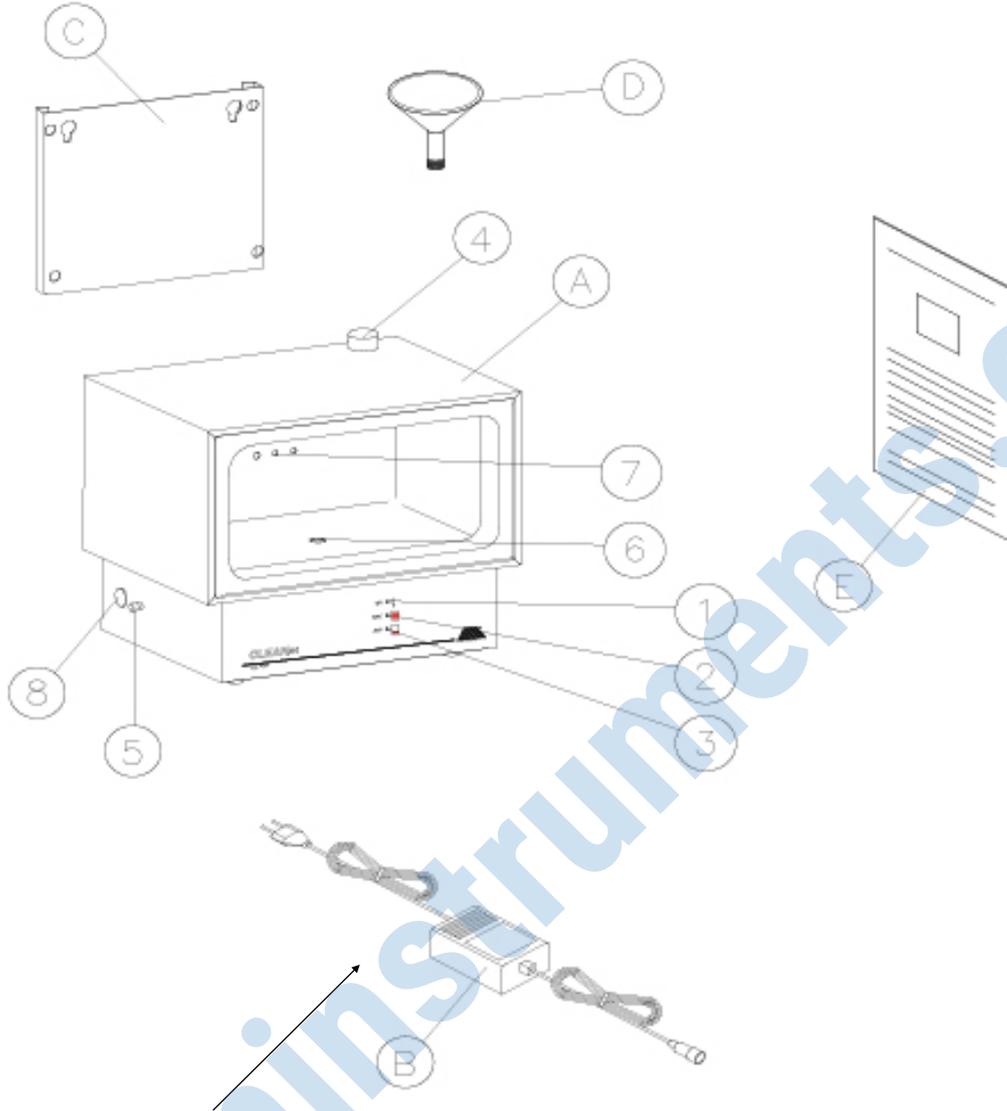non incluso in AMSPC0042
non included in AMSPC0042

<p>AM Instruments srl Via Isonzo, 1/C 20812 Limbiate (MB) Italy</p>	<p>T. +39 02 872892.1 F. +39 02 872892.00 www.aminstruments.com info@aminstruments.com</p>	<p>AM Instruments Tuscany Business Resort Strada di Cerchiaia 28/A 53100 Siena - Italy T. +39 02 872892.1</p>	<p>Sede legale Via Manzoni, 144/E-F 20811 Cesano Maderno (MB) Italy</p>	<p>COMPANY WITH QUALITY SYSTEM CERTIFIED BY DNV GL = ISO 9001:2015 =</p>
--	--	--	--	---

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5.3 - Liquid level indication.

When the liquid level falls below minimum the red light MIN (3) will light up. The CLEANjet can still work normally but each time a hand is introduced, a sound signal will be linked to each spray in order to remind the user to fill up again (see 3.1).

5.4 - Manual running.

The red button (5) can be used for manual running thus excluding the present functions of the optical sensor. Do not operate the CLEANjet in manual mode for a long time to avoid the accumulation and leakage of liquid from the CLEANjet.

In case of liquid leakage, intervene as prescribed on the cleaning solution safety data sheet.

5.5 - Cleaning of CLEANjet.

Disconnect the 230V supply voltage (B), wipe the external and internal surfaces with a cloth soaked with the same cleaning solution used in the CLEANjet.

5.2 - Use of the CLEANjet.

Now the CLEANjet is ready for normal running. Introduce one hand at a time making sure it is in the upper part of the unit and that the fingers are slightly bent to activate the optical sensor (7). The atomization will automatically start up for a preset time (2/3 seconds).

Wait until the end of the cycle and repeat the same operation with the other hand, then rub both hands to distribute the cleaning solution

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7 - WARRANTY

The product and accessories are covered by a guarantee of good functioning for a period of 12 months from the date of sale by AM Instruments (date of the DDT).

The warranty includes the complete and free replacement of the parts deemed, in our opinion, defective in origin, as well as the labor necessary for repair or replacement.

The warranty is only valid for goods delivered ex works.

The costs of transport, those relating to any packaging and charges for services requested by the Customer at their premises (such as tests, checks, inspections for assistance, etc.) are the responsibility of the applicant.

The warranty is not valid if:

- the product and its accessories have not been correctly used, according to the instructions contained in this manual;
- any tampering or modifications have been made by the user or by third parties not authorized by us.

Any damage resulting from lack of or reduced functionality of the appliance, however occurred, does not give the right to any compensation.

To purchase accessories and / or supply spare parts, contact AM Instruments.

In each contact, always indicate the serial number of the appliance and the date of purchase.

AM Instruments srl Via Isonzo, 1/C 20812 Limbiate (MB) Italy	T. +39 02 872892.1 F. +39 02 872892.00 www.aminstruments.com info@aminstruments.com	AM Instruments Tuscany Business Resort Strada di Cerchiaia 28/A 53100 Siena - Italy T. +39 02 872892.1	Sede legale Via Manzoni, 144/E-F 20811 Cesano Maderno (MB) Italy	COMPANY WITH QUALITY SYSTEM CERTIFIED BY DNV GL = ISO 9001:2015 =
--	--	--	--	---