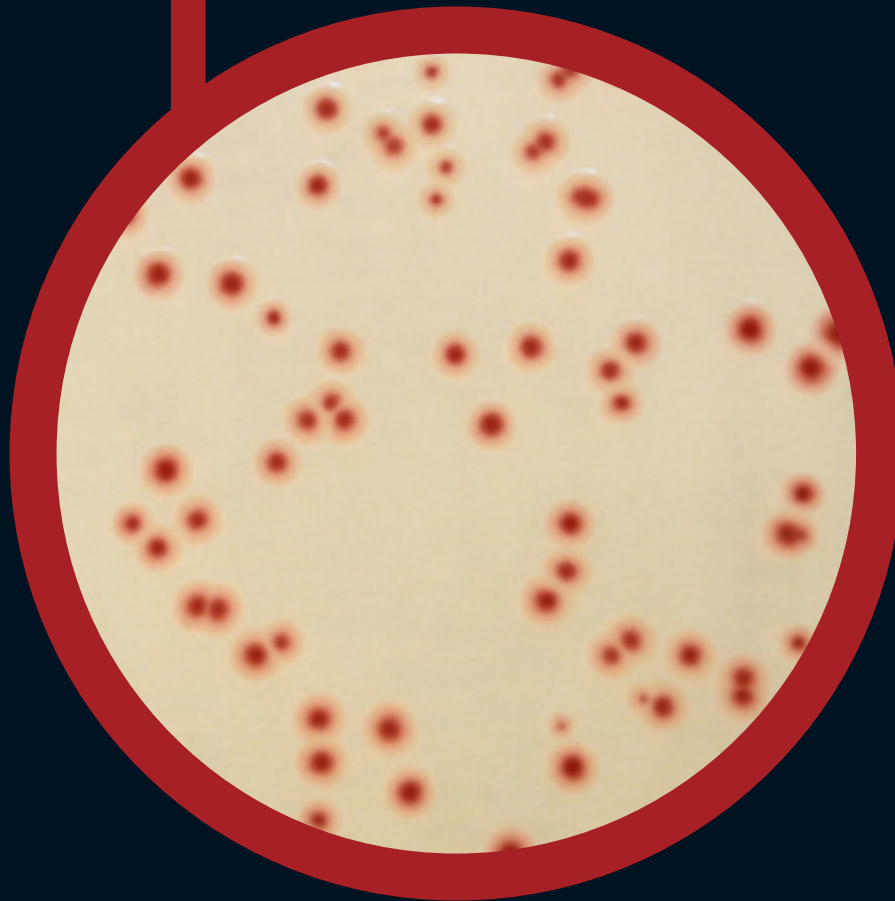
● CHROMagar™
Campylobacter

For detection, differentiation and enumeration
of thermotolerant *Campylobacter*


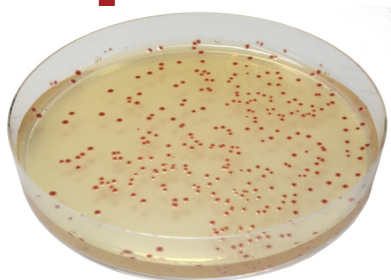
Plate Reading

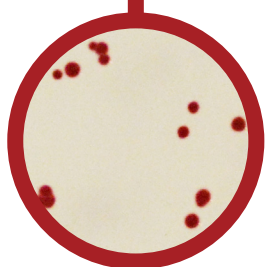- *Campylobacter jejuni, coli, lari*
→ red
- Other bacteria
→ blue or inhibited

For detection, differentiation and enumeration of thermotolerant *Campylobacter*

Background

"*Campylobacter* bacteria are a major cause of foodborne diarrhoeal illness in humans and are the most common bacteria that cause gastroenteritis worldwide. In developed and developing countries, they cause more cases of diarrhoea than foodborne Salmonella. The high incidence of *Campylobacter* diarrhoea, as well as its duration and possible sequelae, makes it highly important from a socio-economic perspective. In developing countries, *Campylobacter* infections in children under the age of two years are especially frequent, sometimes resulting in death." World Health Organisation (WHO) – fact sheet N°255

Campylobacter spp. are fastidious bacteria that may be difficult to recover due to suboptimal specimen transport and/or storage conditions and lack of proper culture procedures. Several culture media formulations have been developed, some blood-based other including charcoal. Each of these media have shown to be a mediocre compromise between specificity and sensitivity.

With CHROMagar™ *Campylobacter*, no compromises! This medium allies the highly specific and easy to read chromogenic technology with an unrivalled growth agar base.

Medium Performance

- 1 IN ACCORDANCE TO ISO 10272-1:2006 REQUIREMENTS**
as a selective medium for the detection of *Campylobacter*
- 2 EASY READING**
36-48h incubation at 42°C.
The intense red colored colonies on a translucent agar facilitates the reading compared to charcoal based agar.
- 3 HIGH RECOVERY RATE ~100%**
Enumerating *Campylobacter* has never been so simple and so reliable.
- 4 HIGHLY SELECTIVE**
In addition, a reinforced selectivity allows for a very clean plate, even with heavily flora loaded samples.

Medium Description

Powder Base	Total	51.2 g/L
	Agar	15.0
+ Supplement (included in the pack)	Peptones and yeast extract	25.0
	NaCl	9.0
	Chromogenic and selective mix	2.2
	Storage at 15/30°C - pH: 7.4 +/-0.2	
	Shelf Life	2 years
	Specific Powder supplement	0.21 g/L
	Storage at 2/8°C	
	Aspect: Powder Form	
	Shelf Life	2 years

Usual Samples	Clinical: Stools Industrial: Food, environmental samples
Procedure	Direct streaking. Incubation 36-48 h at 42°C Micro-aerophilic conditions.
Scientific Publications on this product: available on www.CHROMagar.com Please read carefully the instructions for use (IFU document) available on www.CHROMagar.com	

Quality Control Strains

<i>C.jejuni</i> ATCC® 33291	red
<i>C.coli</i> ATCC® 33559	red
<i>C.lari</i> ATCC® 35221	red
<i>E.faecalis</i> ATCC® 29212	inhibited
<i>C.albicans</i> ATCC® 60193	inhibited
<i>E.coli</i> ATCC® 25922	inhibited

ATCC® is a registered trademark of the American Type Culture Collection

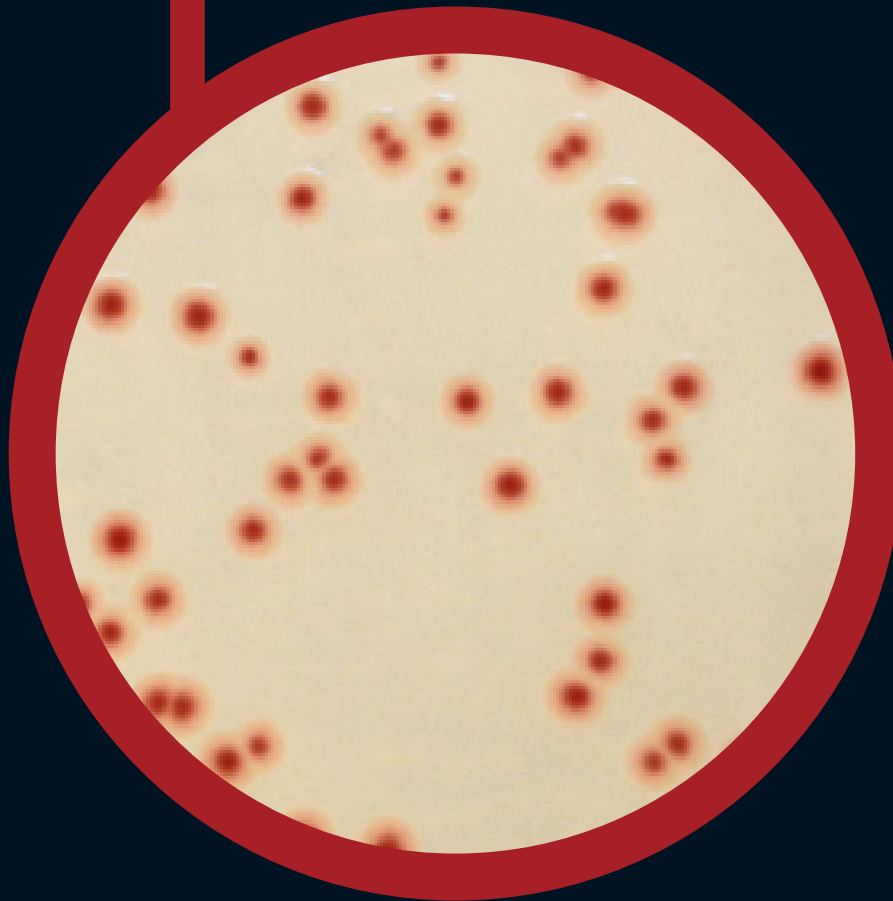
Order References

Please use these product references when contacting your local distributor:

- 5000 ml pack CP572
- 25L pack CP573-25
- (Product = base powder (B) + supplement (S))

Manufacturer: CHROMagar
4 place du 18 juin 1940 75006 Paris - France
Email: CHROMagar@CHROMagar.com
Website: www.CHROMagar.com
Find your nearest distributor on
www.CHROMagar.com/contact

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