

Diagnostics & Ready Prepared Media is our Culture!



# Ready-to-use InstaMedia pouches

for food pathogen testing

We offer a wide range of sterile dehydrated culture media in preweighed and gamma-irradiated bags for direct food pathogen testing. No more media preparation steps, just add sterile water, your food sample and incubate or make your dilutions.



# Ready-to use InstaMedia pouches for food pathogen testing

We offer a wide range of sterile dehydrated culture media in preweighed and gamma-irradiated bags for direct food pathogen testing.

No more media preparation steps, just add sterile water, your food sample and incubate or make your dilutions.

- ▼ Save time and money
- ▼ Save lab space
- ▼ Availability of the shelf
- ▼ Possibility to adapt or customize filling volume based on your handling.

Non-selective enrichment broths available for:

- ▼ Salmonella and Enterobacteriaceae  
*Buffered Peptone Water (conform ISO 21528-1, ISO 6579, ISO 6887-1)*
- ▼ Listeria  
*Demi-Fraser Broth (conform ISO 11290-1)*





# Instructions

1

Take your InstaMedia bag



Weigh your food-sample

2

3

Add food-sample to our pre-weighed bags



Add (pre-warmed) sterile water to the bag

4

5

In case of stomacher bag, pre-dissolve gently and place into stomacher system and incubate.



Visit our website for more information about ready-to-use culture media for secondary enrichments broth.

Product Name	Product Format	Product Code
BPW InstaMediA	Powder for 225 ml, Aluminium bag 700 ml, gamma irradiated	K252I225UR
BPW InstaMediA	Powder for 900 ml, Aluminium bag 1750 ml, gamma irradiated	K252I900JC
BPW InstaMediA	Powder for 2250 ml, Aluminium bag 7500 ml, handle, gamma irradiated	K252I2250JE
BPW InstaMediA	Powder for 2700 ml, Aluminium bag 7500 ml, handle, gamma irradiated	K252I2700JE
BPW InstaMediA	Powder for 3375 ml, Aluminium bag 7500 ml, handle, gamma irradiated	K252I3375JE
BPW InstaMediA	Powder for 3375 ml, 10 L bag, black screw cap, pressure filter, gamma irradiated	K252I3375JJ
BPW InstaMediA	Powder for 6750 ml, 10 L bag, black screw cap, pressure filter, gamma irradiated	K252I6750JJ
Demi-Fraser InstaMediA	Powder for 225 ml, Aluminium bag 700 ml, gamma irradiated	K198I225UR
Demi-Fraser InstaMediA	Powder for 2250 ml, Aluminium bag 7500 ml, handle, gamma irradiated	K198I2250JE
TSB InstaMediA	Vial 20 ml, grey septum/open felscap, gamma irradiated, 100 pcs	K240I010TD
TSB, 4x Concentrated InstaMediA	Vial 20 ml, grey septum/open felscap, gamma irradiated, 100 pcs	K656I010TD
TSB, 10x Concentrated InstaMediA	Vial 20 ml, grey septum/open felscap, gamma irradiated, 100 pcs	K554I010TD
TSB, 2x Concentrated InstaMediA	Vial 50 ml, grey septum/open felscap, gamma irradiated, 50 pcs	K878I050QY

Please contact us if you need a specific product, we are meeting many customers' unique requests.



**bioTRADING Benelux BV**  
 Communicatieweg 7 | 3641 SG Mijdrecht  
 The Netherlands  
 T +31(0)297-286848  
 E sales@biotrading.com

**bioTRADING Belgium BVBA**  
 Haachtsebaan 143 | 3140 Keerbergen  
 Belgium  
 T +32(0)15-500100  
 E info@biotrading.be

Visit our website for more information: [www.biotrading.com](http://www.biotrading.com)