

Fraser Broth Complete

TN 1035 500 g



Intended use

Used for the selective enrichment of *Listeria* spp. from foodstuffs (according to § 64 German Food and Feed Code) and from environmental material.

Form

Powder

Colour

Bright beige

Storage

Dry, tightly closed, at 10...25 °C.

Shelf life

3 years

Typical composition

Component	g/l
Meat peptone	5
Casein peptone	5
Meat extract	5
Yeast extract	5
Sodium chloride	20
Disodium hydrogen phosphate	12
Potassium dihydrogen phosphate	1.35
Esculin	1
Lithium chloride	3
Ammonium iron (III) citrate	0.5
Acriflavin hydrochloride	0.025
Nalidixic acid	0.02

Directions

Dissolve 57.9 g in 1 litre distilled water and autoclave at 121 °C for 15 minutes. Precipitates may be formed after autoclaving. Such precipitate formation has no effect on the test result.

Final pH at 25 °C

7.2 ± 0.2

Microbiological quality control

Test strain	CFU/ 10 ml	Smear (10 µl) Fraser Broth on Palcam Agar after 48 ± 3 h, 36 ± 1 °C
<i>Listeria monocytogenes</i> Serovar 1/2 ATCC 15313	10-100	> 10 colonies per tube
<i>Listeria monocytogenes</i> Serovar 4a ATCC 19114	10-100	> 10 colonies per tube
<i>Listeria monocytogenes</i> Serovar 4b ATCC 19115	10-100	> 10 colonies per tube
<i>Listeria innocua</i> Serovar 6a ATCC 33090	10-100	> 10 colonies per tube
<i>Enterococcus faecalis</i> ATCC 29212	100-1000	no growth

Date of revision: 2014/03/20

Manufacturer: sifin diagnostics gmbh | Berliner Allee 317-321 | 13088 Berlin | Germany | T: +49 30 927 030-0 | F: +49 30 927 030-30 | info@sifin.de | www.sifin.de